
**CELESTRON**  
**MICRO FI**  
Wi-Fi Handheld Microscope  
STREAM • CAPTURE • SHARE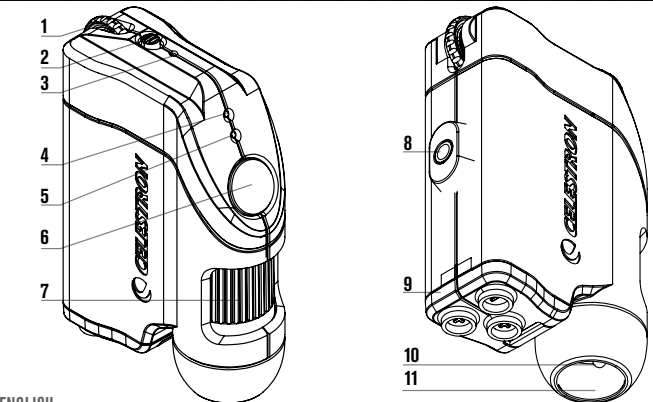

ENGLISH

**I. COMPONENTS**

- 1. ILLUMINATOR ADJUSTMENT KNOB
- 2. ON/OFF SWITCH
- 3. RESET BUTTON
- 4. BATTERY INDICATOR LED
- 5. POWER AND WI-FI SIGNAL INDICATOR LED
- 6. SNAPSHOT BUTTON
- 7. FOCUS WHEEL
- 8. MOUNT FOR TRIPOD OR STAND (1/4-20 STANDARD THREAD)
- 9. BATTERY DOOR
- 10. 6 LED RING ILLUMINATOR
- 11. CLEAR LENS COVER

**II. Getting Started**

- Using your smart device, visit the Apple App Store or Google Play to download the FREE Micro Fi app.
- While the app is downloading, prepare Micro Fi by opening the battery door and inserting 3 AA batteries in the battery compartment (fig. 1).

**NOTE:** Follow the markings on the outside of the battery door to ensure the batteries are inserted correctly (fig. 2).

- Slide the ON/OFF switch (2) to the ON position. The green LED (5) will illuminate.
- Allow 30 seconds for the Wi-Fi network to boot up.

**NOTE:** If Battery indicator LED (4) is illuminated red, the batteries are low on power and will need to be changed out.

**NOTE:** All Micro Fi microscopes create their own Wi-Fi network. Each is named "MICRO-FI" followed by 5 random digits to differentiate the networks. The five digits for your Micro Fi's network are printed on a sticker inside the battery door (fig.3).

- To join the network, enter the default password: 12345678
- When connection to the network is successful, the green Wi-Fi signal LED (5) will blink.
- On your smart device, open the Micro-Fi app.

**NOTE:** You can change the network's name and password in the "Settings" section of the Micro Fi app. To reset name and password back to default, use a paper clip to press and hold the reset button (3) for 3 seconds.

**NOTE:** If the live stream does not start automatically, touch the "ON" button on the screen.

**III. Using the App**

- ICONS
  - **ON/OFF:** Turns live stream on and off.
  - **PAUSE:** Freezes the live feed for shake-free images. (NOTE: This feature available using iOS app only.)
  - **SNAPSHOT:** Takes a still image.

- **RECORD/STOP:** Starts and stops video recording.
- **PLAY:** Displays all saved videos for playback.
- **SETTING:** Allows you to change the Wi-Fi network name and password. On iOS devices, you can also adjust the following image settings: • Brightness • Contrast • Hue • Saturation

**IV. Basics**

- The closer the microscope is to the specimen, the higher the magnification will be.
- To bring a specimen into focus, turn the focus wheel (7) all the way to the right (counter clockwise), then turn it to the left (clockwise) slowly until the specimen comes into sharp focus.
- To obtain the highest magnification of 80x, place the microscope on top of the specimen, with clear lens cover (11) touching it. Continue to turn the focus wheel clockwise until the specimen is in focus.
- Use illuminator adjustment wheel (1) to adjust the light if necessary.
- Use the snapshot button (6) to take a still image. Alternatively, you can press on your app.
- The snapshot images will be saved to your smart device
  - Camera Roll for iPhone, iPad and iPod Touch
  - Gallery for Android devices
- Videos must be taken using the app and pressing .
- Videos are saved to the app and can be viewed by pressing .

**V. Specifications:**

<b>Magnification Ratio</b>	1x to 80x
<b>Digital Camera</b>	.3MP, Color CMOS 1/4"
<b>Frame Rate</b>	15fps (VGA)
<b>Support Video Code Protocol</b>	MPEG4
<b>Shutter trigger</b>	On microscope housing or through app
<b>Lens</b>	2-element IR coated lens
<b>Focuser Range</b>	Manual focus from 10mm to infinity
<b>Illuminator</b>	Upper, 6 bright white adjustable LEDs
<b>White Balance - Auto</b>	<b>Exposure - Auto</b> <b>Gain Control - Auto</b>
<b>Operating Temperature</b>	0° C to 50° C (32° F to 122° F)
<b>Operating Systems Supported</b>	iOS, Android
<b>Support Protocol</b>	2.4GHz Wi-Fi IEEE802.11b/g/n
<b>Security</b>	WPA
<b>Wi-Fi Range</b>	30 feet (without obstructions)
<b>Power Consumption</b>	3.6V/ 440mA
<b>Battery</b>	AA Batteries- QNTY:3 (Not included)
<b>Dimensions</b>	3.15" x 1.36" x 4.39" (80.10mm x 34.50mm x 111.6mm)
<b>Weight</b>	121g/42. Oz.

**Warranty:** Your microscope has a two year limited warranty. Please see Celestron's website for detailed information on all Celestron microscopes at [www.celestron.com](http://www.celestron.com).

**FCC Note:** This equipment has been tested and found to comply with the limits for a Class B digital device, pursuant to part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference in a residential installation. This equipment generates, uses, and can radiate radio frequency energy and, if not installed and used in accordance with the instructions, may cause harmful interference to radio communications. However, there is no guarantee that interference will not occur in a particular installation. If this equipment does cause harmful interference to radio or television reception, which can be determined by turning the equipment off and on, the user is encouraged to try to correct the interference by one or more of the following measures:

- Reorient or relocate the receiving antenna.
- Increase the separation between the equipment and receiver.
- Connect the equipment into an outlet on a circuit different from that to which the receiver is connected.
- Consult the dealer or an experienced radio/TV technician for help.

This device complies with Part 15 of the FCC Rules. Operation is subject to the following two conditions: (1) this device may not cause harmful interference, and (2) this device must accept any interference received, including interference that may cause undesired operation.

**RF WARNING**

This equipment complies with FCC RF exposure limits set forth for an uncontrolled environment. This equipment must be installed and operated with a separation distance of at least 20 cm from all persons. Apple, the Apple logo, iPhone, iPod touch and iPad are trademarks of Apple Inc., registered in the U.S. and other countries. App Store is a service mark of Apple Inc.

**This product is designed and intended for use by those 14 years of age and older. Product design and specifications are subject to change without prior notification.**

**FOR COMPLETE SPECIFICATIONS AND PRODUCT INFORMATION, VISIT: WWW.CELESTRON.COM**  
Celestron, LLC. • Torrance, CA 90503 U.S.A. Tel: 800.421.9649 ©2014 Celestron • All rights reserved. • Printed in Taiwan

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