## DIGITAL MICROSCOPE IMAGER

INSTRUCTION MANUAL  
MODEL #44421

ENGLISH

## ***INTRODUCTION***

Congratulations on your purchase of the Celestron Digital Microscope Imager. Your Imager is a precision optical instrument made of high quality materials to ensure durability and long life, and designed to give you a lifetime of pleasure with minimal maintenance. It is miniature in size – 2.5" (64 mm) long x 0.9" (22 mm) in diameter and weighs just 1.5 oz (42 g) but powerful in its application with a 2 MP sensor for the digital camera.

This Imager can be used for virtually anything that traditional microscopes are used for: Hobbyists, Educators, Medical Labs, Industrial Inspection, Engineering Applications, Teachers, Students, Science Applications, Doctor's Offices, Police Agencies, Government Testing, and general use by consumers – except you can view the images on a computer screen and take snapshots or videos of the images.

The Digital Microscope Imager will work with virtually any microscope (biological and stereo) as long as the eyepiece barrels are of the universal sizes of 23 mm and 30 mm or similar. You replace the eyepiece in the microscope with the Imager.

Before attempting to use your Imager, please read through the instructions to familiarize yourself with the functions and operations to make usage easier for you.

With the included software, you can observe the magnified images or capture video or take snapshots with Microsoft Windows XP, Vista, and 7 operating systems. You can also use the microscope with other image capture software that you may have on your computer. If you are using a MAC, you can take videos and snapshots but you need image/photo capture software for MAC. You will also need a CD/DVD drive as well as an open USB port.

## Setting Up Your Digital Microscope Imager

1. Carefully remove the Imager, Adapter, and CD-ROM (Digital Microscope Suite) from the box and set them on a table, desk, or other flat surface. The USB cable is permanently attached to the Imager.
2. Remove the plastic bag covering the Imager.
3. The Adapter (image below) will attach to the Imager with three thumbscrews. First back up the threads from the Adapter so they do not protrude through the barrel so there is no interference when inserting the Imager and then insert it.
  - A. The position of the Adapter on the Imager is not critical but the top of the adapter (thumbscrew end) should be about 40 mm (1.6") from the sensor end for optimal performance on most microscopes. Tighten the screws.

4. Remove the eyepiece (just one eyepiece if your microscope uses two eyepieces as with a stereo microscope or binocular head with a biological microscope).
5. Put the Imager into the eyepiece barrel. It is held in place by its own weight and this is ok. However, when moving your microscope with the Imager in the eyepiece barrel, do not turn the microscope sideways or upside down or the Imager will fall out.

## Using the Imager

Before using your Imager for viewing on your computer screen or imaging, you will have to ensure the driver is recognized by your computer. Your Imager is a "Plug and Play" type where the computer should automatically recognize the driver type on Microsoft Windows XP, Vista and 7 Operating Systems as well as Apple MAC version 10.4.9 and later Operating Systems.

Plug the USB cable of the Imager into the USB port of your computer. Each computer system will have slightly different messages. For example, on some Vista OSs when plugged in the message may say "Installing Device Driver Software" and shortly thereafter you may see a message "Your Devices Are Ready to Use". Just follow any instructions shown on the screen.

You will need to install the Digital Microscope Suite software (on Windows OS) to actually view and/or image and prior to doing this you will need to unplug the USB cable until the installation is complete. You can also use other image/photo capture software that you may have on your Windows system if you desire. To use the Imager with MAC operating systems you will need to have image/photo capture software on your computer (like a combination of "iChat" with "Photo Booth", etc.).

The approximate digital power is 100x (on a monitor of 14" (35 mm) to 30" (75 mm) using the Imager. The actual power is determined by the digital power using the imager (100x) times the objective lens power. As an example, if the objective lens being used is 4x, then the power you will be capturing for snapshots or video will be 400x. If using a 40x objective, then the power will be 4000x, etc.

## **Adjusting Light Intensity**

Whether using mirror illumination or electrical illumination, you will have to experiment with how much light will give you the best image. You can also experiment with condensers and diaphragms (if your microscope has these) for optimum adjustment for obtaining the best image.

## **Focusing**

You will focus your microscope with the Imager attached the same as you would for observing. For the best images, make sure the specimen or object is sharply focused. When changing powers, you will have to refocus.

## **Rotating the Imager**

When you are looking at the image on your computer screen, you can rotate the position 360° by rotating the Imager to the position you prefer.

## **Installing the Digital Microscope Suite (DMS) Software**

Your Imager comes with the Digital Microscope Suite CD-ROM which allows you to use it for visual purposes and for imaging with your computer when using Microsoft Windows OSs.

Before installing the software, remove the USB cable of the Digital Camera from your computer and do not connect it until the software installation is complete. Insert the CD-ROM into your CD-ROM drive and it should automatically begin. Messages will appear (each operating system may show something somewhat different) to guide you through the installation process.

If the software did not begin automatically, double click on "(My) Computer" and right click the CD-ROM icon and it should start.

- You will see "Files Currently on the CD" and notice "Digital Microscope Suite" as a folder. Double click on it and then among the choices double click on "Setup" (InstallShield Setup Launcher).

- Then, the InstallShield Wizard will appear. Then select the default choices, “Next”, etc. until “Finish” is seen and click on it.

Your software was properly installed. You will have a desktop icon called “Digital Microscope Suite” but if not that will be the program you select from your installed programs when you want to use the digital microscope Imager.

## Using the Digital Microscope Suite (DMS) Software

Plug in the USB cable of the Imager to your computer and then start DMS. The DMS contains basic image capture and organization for your photos (snapshots) and videos. Click on the DMS icon to begin.

The software is very intuitive and easy to use. You can change settings by using File and Settings – the Video Device should have as a default “Celestron Microscope Imager” but if not please change it so the DMS will function properly.

You can take snapshot images by using the snapshot button on the microscope (only with Windows OS) or you can also use the keyboard by clicking enter on “Capture an Image”. For videos you need to start/stop on the icon shown on the screen.

**You are now ready to begin your digital adventure!**

## **CARE, MAINTENANCE, AND WARRANTY**

Your Imager is a precision optical instrument and should be treated with care at all times. Follow these care and maintenance suggestions and your Imager will need very little maintenance throughout its lifetime.

- Store the Imager in a dry and clean place.
- Be very careful if using your microscope in direct sun light to prevent damage to the Imager or your eyes.
- Never point the Imager towards the sun or the sensor can be damaged and cease working.
- Clean the outside surfaces with a moist cloth.
- Never clean optical surfaces with cloth or paper towels as they can scratch optical surfaces easily.
- Blow off dust with a camel's hair brush or air blower off the optical surfaces.
- To clean fingerprints off of optical surfaces, use a lens cleaning agent and lens tissue available at most photo outlets and when cleaning do not rub in circles as this may cause streaks or scratches to occur.
- Never disassemble or clean internal optical surfaces. This should be done by qualified technicians at the factory or other authorized repair facilities.
- Do not disassemble or take apart the Imager or damage can occur.

## **WARRANTY**

Your Imager has a two year limited warranty. Please see the Celestron website for detailed information at [www.celestron.com](http://www.celestron.com).

www.celestron.com

2835 Columbia Street • Torrance, CA 90503 U.S.A.

Telephone: 310.328.9560 • Fax: 310.212.5835

©2012 Celestron

All rights reserved. • Printed in China • 04-12

FCC Note: This equipment has been tested and found to comply with the limits for a Class B digital device, pursuant to part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference in a residential installation. This equipment generates, uses, and can radiate radio frequency energy and, if not installed and used in accordance with the instructions, may cause harmful interference to radio communications. However, there is no guarantee that interference will not occur in a particular installation. If this equipment does cause harmful interference to radio or television reception, which can be determined by turning the equipment off and on, the user is encouraged to try to correct the interference by one or more of the following measures:

- Reorient or relocate the receiving antenna.
- Increase the separation between the equipment and receiver.
- Connect the equipment into an outlet on a circuit different from that to which the receiver is connected.
- Consult the dealer or an experienced radio/TV technician for help.

This product is designed and intended for use by those 14 years of age and older.

Product design and specifications are subject to change without prior notification.

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