DIGITAL **MICROSCOPE IMAGER**

SOFTWARE MANUAL

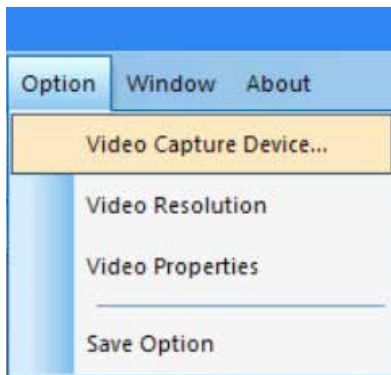
MODEL #44423

Note: The contents of this manual and updates are subject to change without notice.


1. OPERATING INSTRUCTIONS

A. SOFTWARE PRE-SETS


1. When running the S-Viewer software for the first time, taking the time to make a few pre-set choices will make future use easier and more effective.
2. When more than one video device is available on the computer, the software allows you to select the target video device you wish to use.
3. Select **Video Capture Device** from the **Option** tab in the main toolbar.

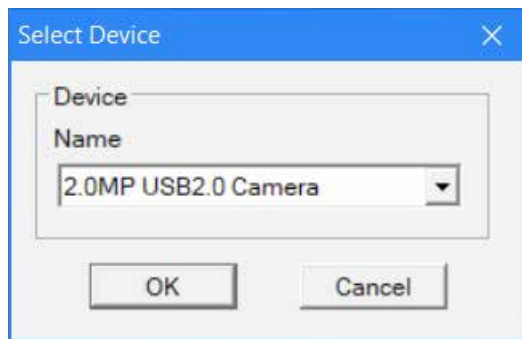
4. When the dialog box **Select Device** appears, you can select the target video device in the list of devices.

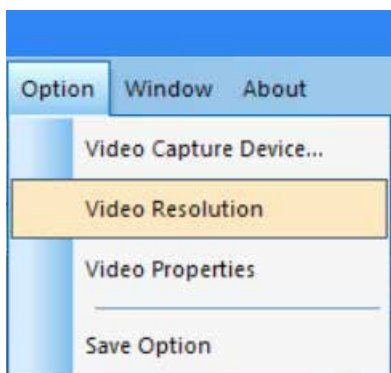5. Next, set the video resolution. To begin, select **Option**→**Video Resolution**.

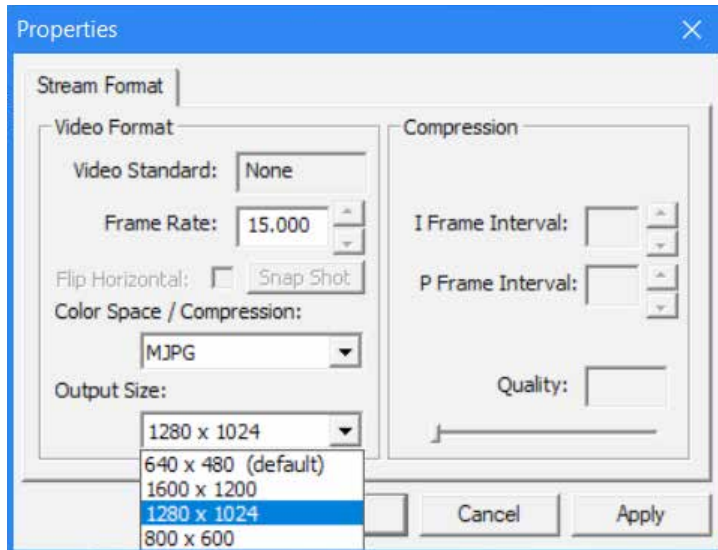
6. When the dialog box **Properties** appears, you can change the video resolution by choosing from the list of options.


7. Next, set the desired video attributes.

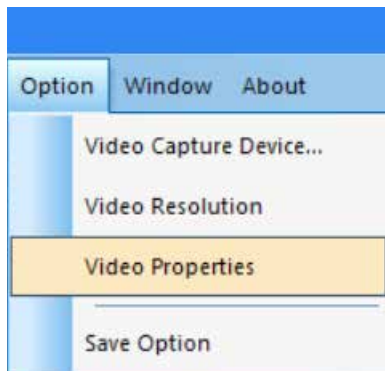8. Select **Option**→**Video Properties** to adjust the basic video capture properties.


9. The **Cap Properties** window will appear.

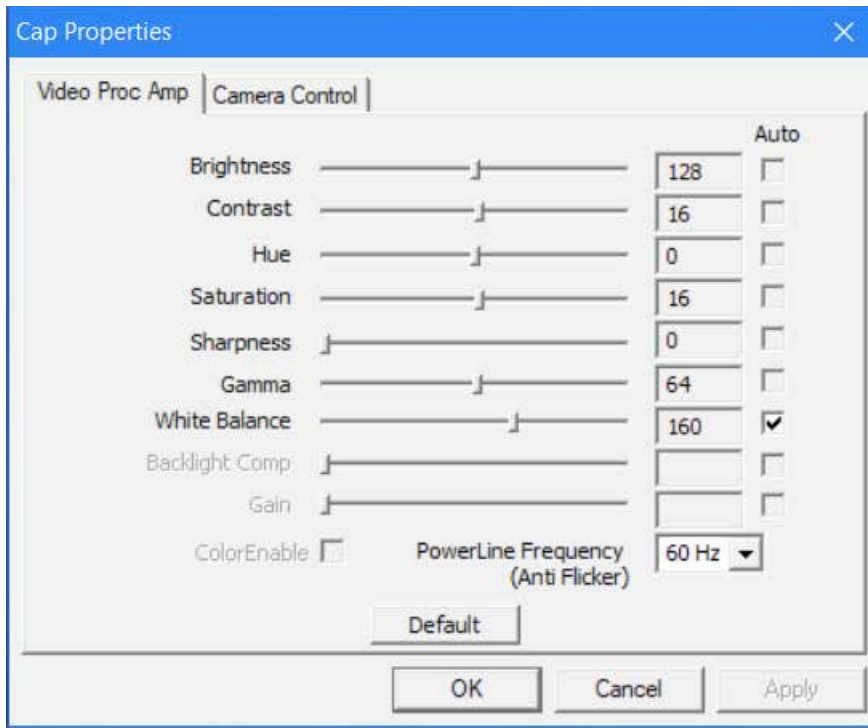
10. The **Video Proc Amp** tab allows you to adjust the video Brightness, Contrast, Hue, and other attributes.

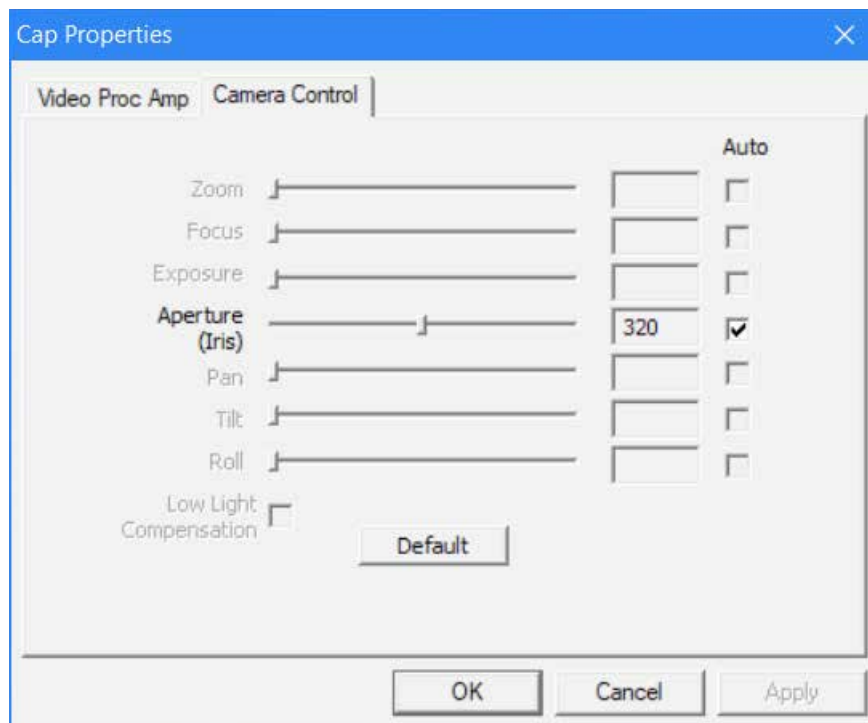
11. The **Camera Control** tab allows you to pre-set the Aperture (Iris) of the camera.

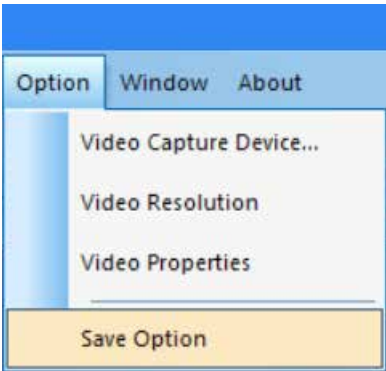
- 12. Finally, select how you would like to save your recordings.
- 13. Select **Option**→**Save Option**, and the Save Option dialog box will pop up.

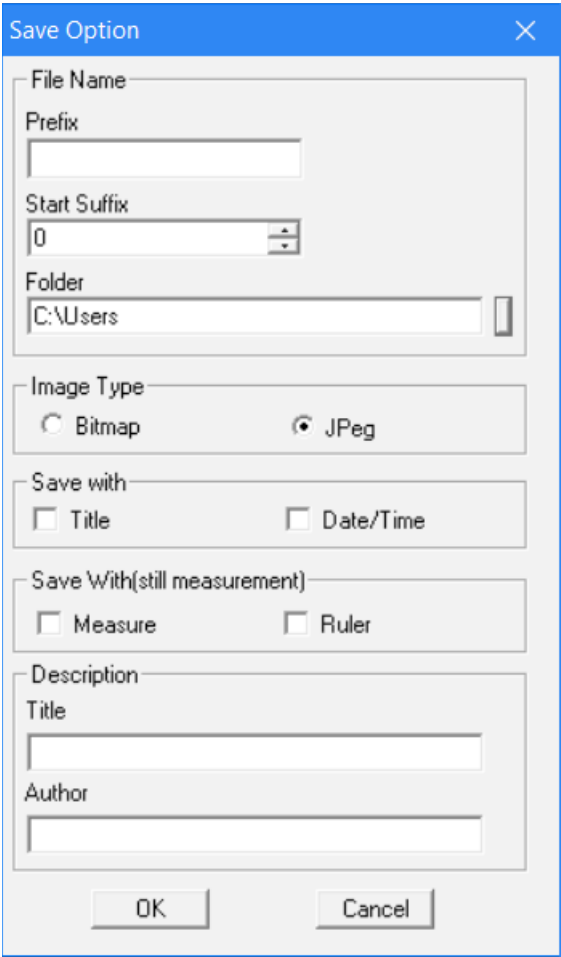- 14. From the **Save Option** window, you can set where and how you would like the recordings to be saved to your computer.

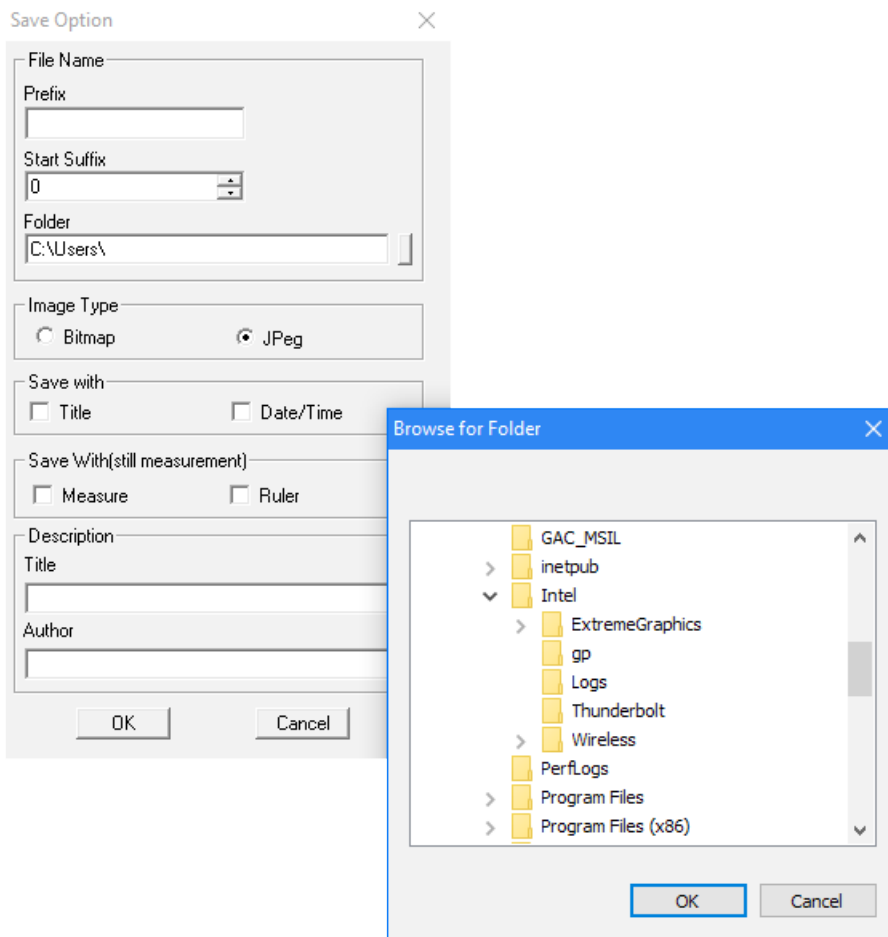15. **Shortcut:** Clicking on the vertical rectangle next to the Folder address bar will bring up the **Browse for Folder** window.


16. Once you have made your selections here, your recordings will always be saved accordingly when you select **File→Save**.

17. **Helpful Tip:** If you ever would like to save a recording to a different location, simply select **File→Save As** and the **Save As** window will appear to enable you to save the file somewhere else on your computer.


B. OPERATION- MENU CONTROLS

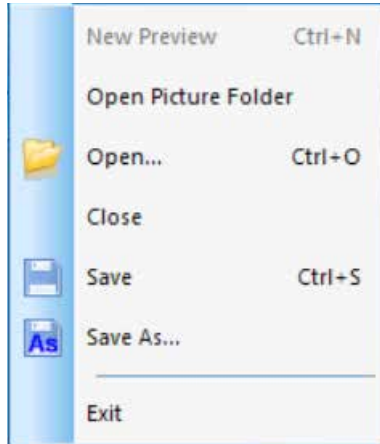FILE MENU


1. **Open Picture Folder:** Opens the folder set as the default location for saving recordings made with the software.
2. **Open:** Allows you to open files from their saved locations.
3. **Close:** Closes the opened videos or images.
4. **Save:** Saves the current video or still image to the default file location. For more information on defining the file location, see **Option→Save Option**.
5. **Save As:** Opens the **Save As** window to enable you to save a recording to a location other than that specified as the default location.
6. **Exit:** Exits the application.

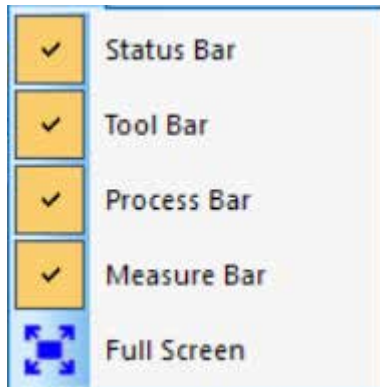VIEW MENU

1. **Status Bar:** Displays or conceals the status bar.
2. **Tool Bar:** Displays or conceals the consequently-used tool bar.
3. **Process Bar:** Displays or conceals the process bar.
4. **Measure Bar:** Displays or conceals the measuring tool bar.
5. **Full Screen:** Displays the video in full screen. (Use the **Esc** key to return to normal view.)

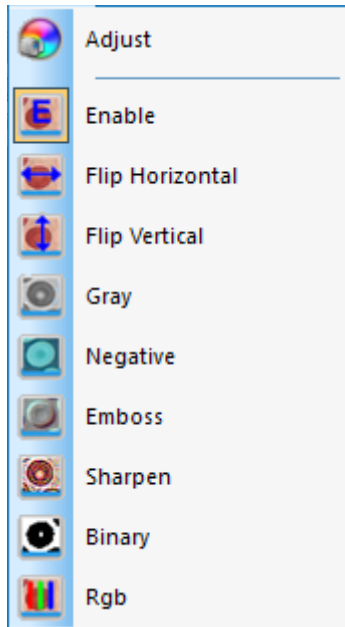
VIDEO MENU


1. **Adjust:** Opens the **Cap Properties** window to allow adjustment of video parameters (Color, Exposure, WB, Frame Rate, Power Frequency, etc.).
2. **Enable:** Activates or deactivates the real-time processing effects (described below) and/or real-time measurement.
3. The rest of the commands in the Video menu are also accessible via the toolbar.

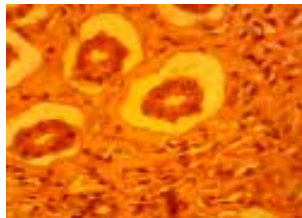The effects of these commands are as follows:


Original Image

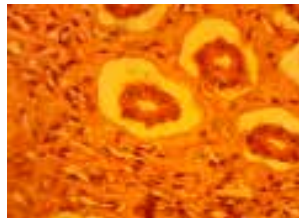
A. Flip Horizontal

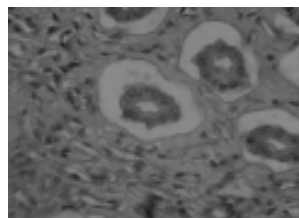B. Flip Vertical

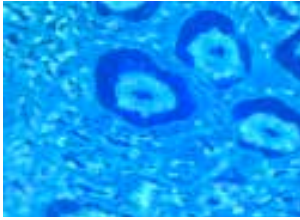C. Gray


D. Negative

E. Emboss

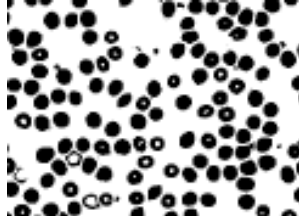
F. Sharpen

G. Binary

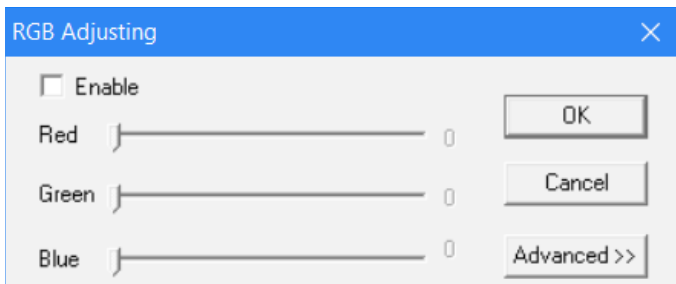
A note on **Binary Video**: You can adjust the threshold value of binary (1 to 255) to display dynamically; simply check the Enable box in the **Binary** window to use or deactivate the binary real-time processing effect.

4. **RGB**: Opens and closes the RGB real-time video control window.


A. To use these RGB controls, simply check the Enable box in the **RGB Adjusting** window.

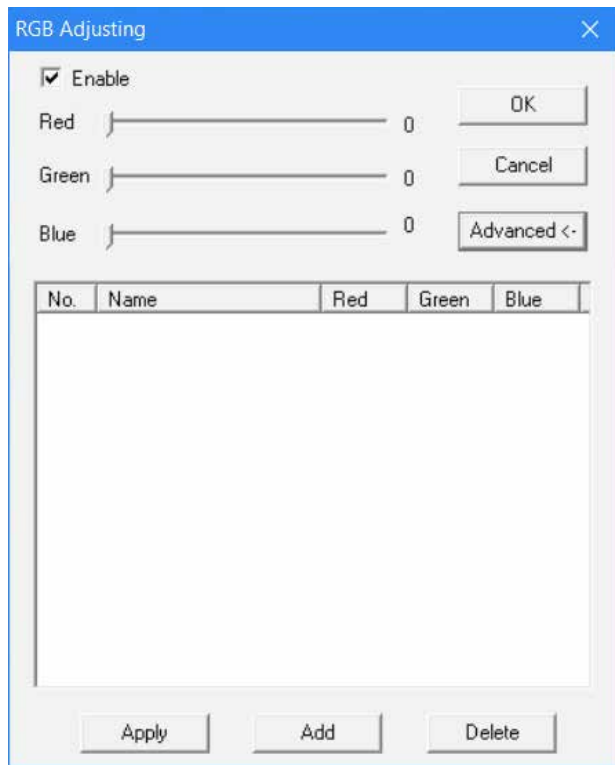B. A control window for advanced RGB adjustments can be accessed through the **Advanced** button.

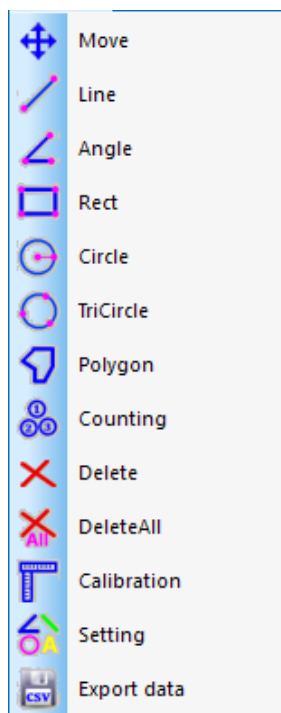- C. Simply check the **Enable** box in this window to activate these controls.
- D. Adjust the red-green-blue slide bars for video tri-color matching.
- E. **Apply**: Selects options from list box and then applies them to the current video.
- F. **Add**: Adds the adjusted red-green-blue value to the database and displays this formula in the list.
- G. **Delete**: Deletes options from the list or database.

MEASURE MENU

1. The commands in the Measure menu can also be accessed in the toolbar under Measure.
 - A. This toolbar is initially positioned vertically and to the right of the image area in the software, but it can be dragged to a horizontal position along the top if desired.


2. The controls in this menu are as follows:

OPERATION TYPE	OPERATION INSTRUCTION	DISPLAYING VALUE
 Move Movement	Selects image to be moved. The image will display in a dotted line. Click and drag to the desired new location.	
 Line Measure point to point	Select the starting point. Click the left button of the mouse and drag to your desired end point.	The length of the line
 Angle Measure an angle	Click three points in different positions with the mouse to form an angle.	The angle's measurement
 Rect Measure a rectangle	Select the top left corner of a rectangle. Click the left button of the mouse and drag to the lower right corner of the rectangle.	The area of the rectangle
 Circle Measure a circle	Select the center of a circle. Click the left button of the mouse and drag to the point on the arc.	The radius of the circle
 TriCircle Measure a circle by three points	Click three points in different positions with the mouse to form a circle.	The radius of the circle
 Polygon Measure the area of a polygon	Select the first point by clicking the left mouse button. Release the button and move to the second point in the polygon. Click the left button to mark that point. Repeat until all points are marked. To close the polygon, move the mouse back to the first point and click the right mouse button.	The area of the polygon
 Counting Identify locations with circled numeral	Select the first point and click the left mouse button. Continue in sequence with all points where markers are desired.	Numbered circles
 Delete Delete an item	Select the item desired to be deleted by clicking the left mouse button.	
 DeleteAll Delete all measurements	Select this command to delete all measurements made.	


3. Special commands:

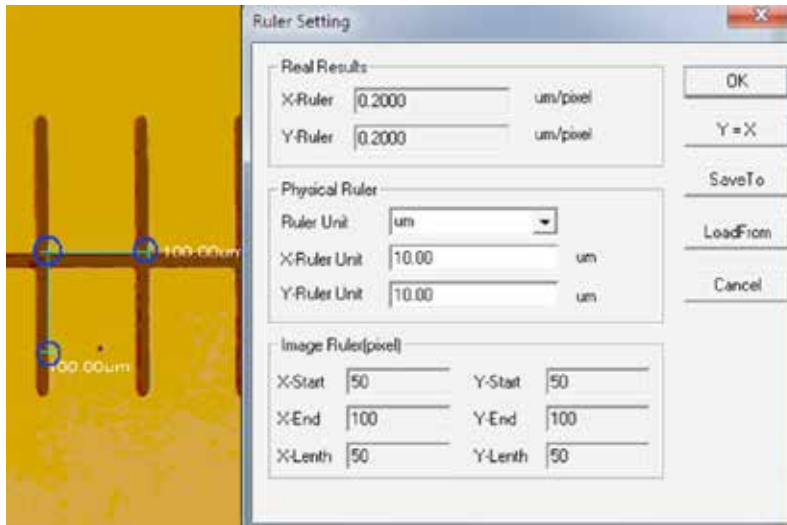- A.  **Calibration** : Selecting the **Calibration** command opens the **Ruler Setting** window. With this window, you can use a standard microscope calibration ruler or other ruler with sufficiently small graduations to calibrate your software's measurement functions.


- i. **Delete Available Units of Measurement:** um/mm/cm/inch.
- ii. **X Direction:** The length along the X axis.
- iii. **Y Direction:** The length along the Y axis.
- iv. **Demarcation Adjustment:** Move the movable points (in the blue circle) of the ruler seen in the image, and the value of trim size (unit length per pixel) will change.
- v. **Y=X:** Makes the X-axis ruler equal to the Y-axis ruler.
- vi. **Save To:** Save the currently demarcated parameter to a file.
- vii. **Load From:** Load a previously demarcated parameter file.

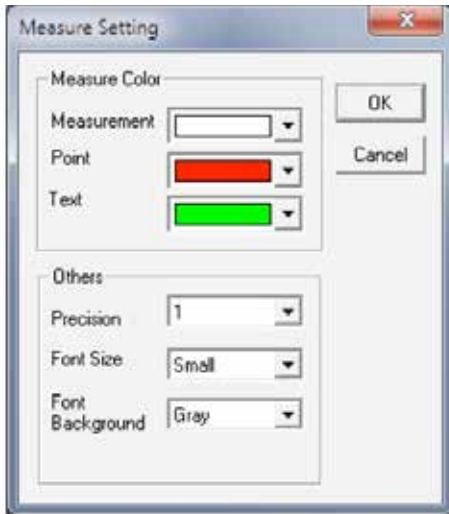
- B.  **Setting** : Selecting the **Measure Setting** command opens a window to enable you to set the colors of measuring lines, movable points, and words.

The image below shows an example of the results of the settings shown in the Measure Setting window illustrated just above:

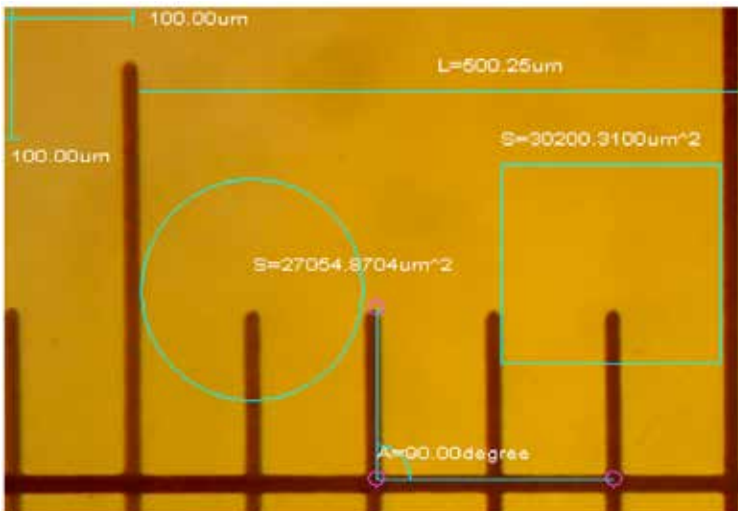
ANNOTATION MENU

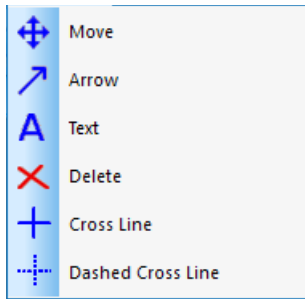- 1. Move:** Selects the added annotation to be moved. The annotation will display in a dotted line. Click and drag to the desired new location.
- 2. Arrow:** Draws an arrow. Click the left mouse button on the location to which the arrow is desired to point then hold and drag in the direction from which you'd like the arrow to originate.
- 3. Text:** Inserts a textbox.

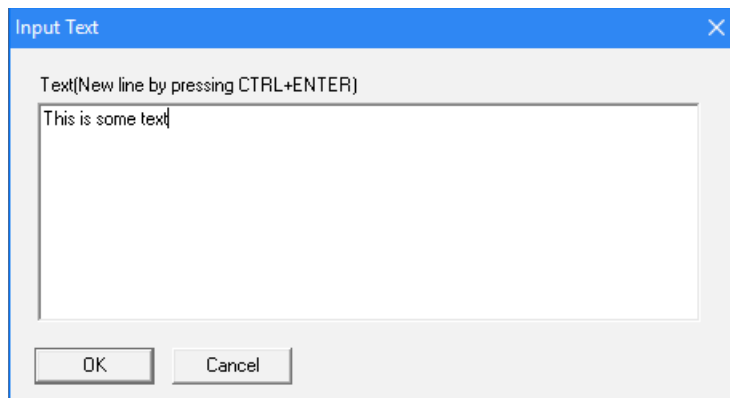- A.** Select the **Text** command.
- B.** The **Input Text** window will open.

- C.** Type your desired text in the box and select **OK**.
 - D.** Your text will be added to the image.
- 4. Delete:** Deletes an annotation.
 - 5. Cross Line:** Overlays a solid X,Y line across the image area.
 - 6. Dashed Cross Line:** Overlays a dashed X,Y line across the image area.
 - 7.** The rest of the commands in the Video menu are also accessible via the toolbar.

STANDARD TOOLBAR

1.  **Open, Save, and Save As:** Same commands as found in the **File** menu.

2.  **Preview:** Activates or returns to live image viewing.

3.  **Pause:** Pauses live image viewing.

4.  **Capture:** Record video images to display in a newly-opened window.

5.  **Interval Capture:** Opens the **Interval Capture** window.

A. **Wait (sec):** Seconds after which the image capture begins.

B. **Interval (sec):** Duration between image captures.

C. **Images:** Number of images to be recorded.

6.  **Record:** Opens the **Record** window.

A. **Path:** Selects the address for saving video files.

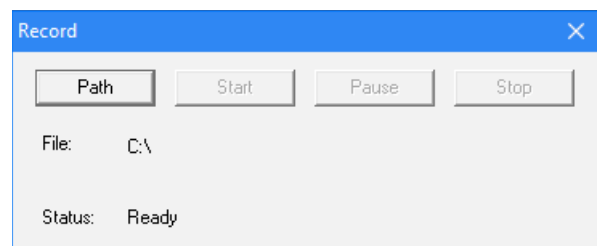
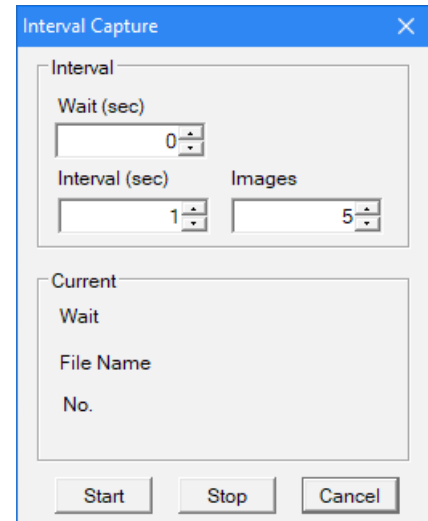
B. **Start:** Start video recording.

C. **Pause:** Pause the video.

D. **Stop:** Stop recording.

7.  **Adjusting Video:** Opens the **Cap Properties** window (the same menu as used to set the video properties in the Software Pre-sets section of this manual).

8.  **Toggle Full Screen:** Displays the image area at full screen. (Use the **Esc** key to return to normal viewing.)

2. TROUBLE-SHOOTING

A. ISSUE: NO PREVIEW (BLANK SCREEN)

Follow these steps:

1. Check to ensure the microscope's illumination source is activated or its mirror properly positioned.
2. Check to ensure that the video device is connected, and the video device driver is installed in Device Manager.
3. Check to ensure the Exposure (Iris) value under **Cap Properties** is normal.
4. Check the dialog box **Select Video Device** to make sure the correct target video device is selected.
5. Close the software and unplug the imager. Then, plug in the imager again and restart both computer and software.

If the unit and software still don't function properly after you have followed all these steps, please contact Customer Service:

Celestron
Customer Service Department
2835 Columbia Street
Torrance, CA 90503

Tel. (310) 328-9560
Fax. (310) 212-5835
Monday-Friday 8AM-4PM PST

3. CARE AND MAINTENANCE

Your Celestron Digital Microscope Imager is a precision optical instrument and should be treated with care at all times. Follow these suggestions and your microscope will need very little maintenance throughout its lifetime.

- + Store your imager in a clean, dry place.
- + Never use your imager where operation of such devices is restricted. Improper use creates the risk of serious accident.
- + Only use your imager within the 23° to 120°F (-5° to 50° C) temperature range. Try to limit sudden temperature changes as this may create moisture in the imager housing.
- + Do not try to access the internal components of your imager or attempt to modify it in any way. Only authorized technicians should perform maintenance or repairs.
- + Keep your imager away from water and other liquids. Never use it in rain or snow. Moisture creates the danger of fire and electric shock.
- + If the exterior of your imager needs cleaning, wipe it with a moist cloth.

4. WARRANTY

Your Celestron Digital Microscope Imager has a two year limited warranty. Please visit the Celestron website for detailed information on all Celestron products: www.celestron.com/pages/warranty.

©2018 Celestron • All rights reserved • celestron.com
2835 Columbia Street • Torrance, CA 90503 U.S.A.
Telephone: 1(800) 421-9649 • Printed in China

FCC Note: This equipment has been tested and found to comply with the limits for a Class B digital device, pursuant to part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference in a residential installation. This equipment generates, uses, and can radiate radio frequency energy and, if not installed and used in accordance with the instructions, may cause harmful interference to radio communications. However, there is no guarantee that interference will not occur in a particular installation. If this equipment does cause harmful interference to radio or television reception, which can be determined by turning the equipment off and on, the user is encouraged to try to correct the interference by one or more of the following measures:

- + Reorient or relocate the receiving antenna.
- + Increase the separation between the equipment and receiver.
- + Connect the equipment into an outlet on a circuit different from that to which the receiver is connected.
- + Consult the dealer or an experienced radio/TV technician for help.

This product is designed and intended for use by those 14 years of age and older. Product design and specifications are subject to change without prior notification.

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