Display:	¾" LCD
Range:	Counts down from 20 hrs.

Resolution: 1 min.

Accuracy: 0.01%

Case:

**SPECIFICATIONS** 

ABS plastic One 3-year AAA battery

Power: Accessories

Supplied: 3-year AAA battery, magnet/clip, Traceable® certificate, instructions

QC Timer Operation 1. Turn the timer on by inserting the battery. To insert the battery, slide the battery cover unit. Place the supplied battery inside. Replace the battery cover.

2. To set the desired time, press the HR and MIN buttons. Each press advances the time 1 hour or 1 minute, respectively. To rapidly advance

located on the back of the unit away from the

counting down. The blinking colon (:) indicates the time is running. 4. To temporarily stop the countdown time, press

the time, press and hold the buttons.

the START/STOP button once. To resume countdown, press the START/STOP button again.

3. Press the START/STOP button once to begin

**Battery Information** 

CLEAR.

Please replace the battery in the case of any operational difficulties or lack of proper functionality for any reason. Slide the battery cover located on

When the timer reaches zero, an alarm will

button to silence the sounding alarm.

6. When the timer is counting down, clear the

sound for 1 minute. Press the ALARM STOP

display by pressing the START/STOP button

to stop the countdown. Then press the CLEAR

button. If the timer is not counting down, press

the START/STOP button *twice* and then press

the old battery and replace it with a new AAA battery. Replace the battery cover. Incorrectly installed batteries may damage electronics.

the back of the unit away from the unit. Remove

Replacing the battery with a new fresh battery will solve most difficulties.

TRACEABLE® QC TIMER INSTRUCTIONS